

InstiSHAKE™ HEK Supplement

Product Code: RS-2325

Description


InstiSHAKE™ HEK is a chemically defined, animal component free, cell culture supplement designed to support the growth of suspension HEK cells in shaking culture. Specifically, InstiSHAKE HEK supports HEK cells during transition from static to shaking culture conditions. InstiSHAKE HEK is supplied as a 20x concentrated solution which should be diluted to a final concentration of 1x in cell culture medium.

Intended Use

FOR RESEARCH USE ONLY. NOT INTENDED FOR HUMAN, ANIMAL, DIAGNOSTIC, OR THERAPEUTIC USES.

Storage Conditions

Store at 2-8°C. Do not freeze.

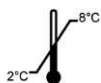

Shelf-Life Information

InstiSHAKE HEK is stable for 18 months from the date of manufacture (see Certificate of Analysis on our website - <https://www.aicompanies.com/support/certificates-of-analysis/>).

Directions for Use

1. Gently invert the InstiSHAKE HEK vial several times (do not shake).
2. Prepare base cell culture media in accordance with in-house practices.
3. Add InstiSHAKE HEK to the cell culture medium to give a final concentration of 1x.

Example: Add 25ml InstiSHAKE HEK to 475ml of cell culture medium.